WILD
HEERBRUGG

SERVICE MANUAL

M420

PHOTOMAKROSKOP

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These instructions are intended for service personnel trained by Wild only. They comprise service level 3 only.

Those using these instructions are required to be familiar with the specific instrument, in addition to having an adequate basic knowledge of mechanics and geometrical optics, together with precision-engineering skills.

- Service level 1: Functional checks, putting into service, maintenance and adjustments in accordance with operating manual
- Service level 2: More extensive maintenance work such as cleaning, lubrication readjusting; simple servicing
A workshop with the necessary special tools is required
This service level also includes service level 1
- Service level 3: All types of repair work
A workshop fully equipped with a complete set of tools is required
This service level also includes service levels 1 and 2

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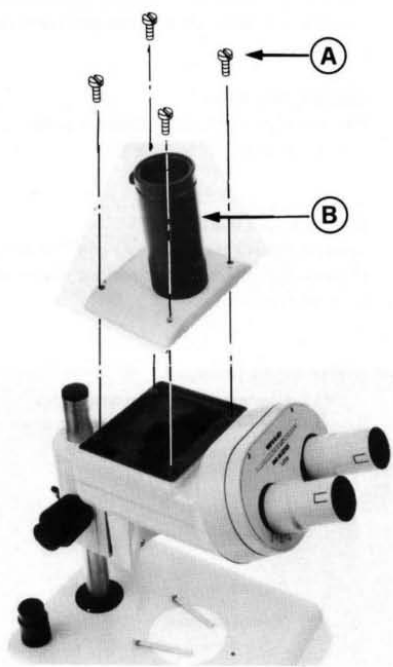
Caution:

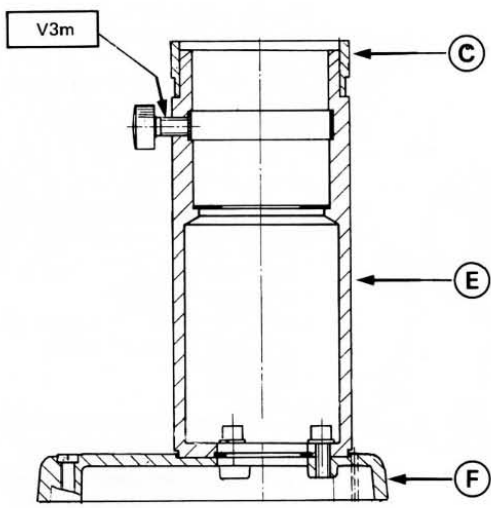
The photo tube (B) serves as a reference for the adjustment. When separating the carrier tube (E) from cover (F), or removing adjusting ring (C), the photo tube must be adjusted immediately before continuing the disassembly. See pages 18,19.

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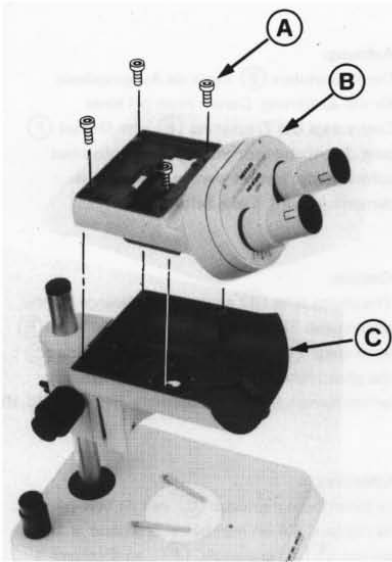
Assembly instruction:

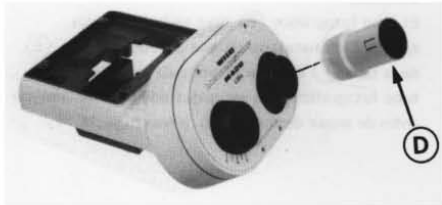
Press upper part (B) against reference surface (C) and fit it symmetrical to lower part.

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Assembly instruction:

- Cover disk (G) must be on top of separating metal-sheet (H)
- For adjusting the eyepiece (D) tubes see adjustment instructions

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
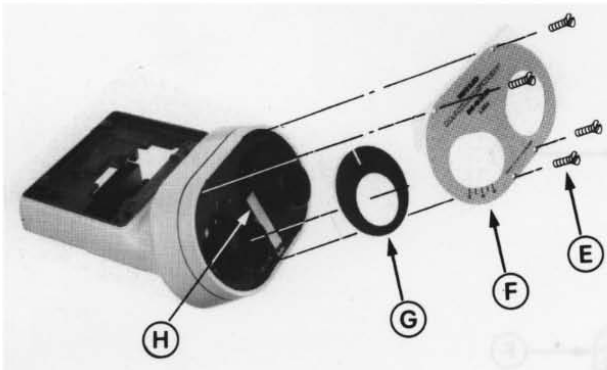
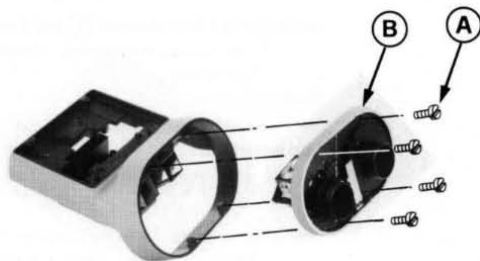
Assembly and adjustment of front plate (B) see adjustment instructions

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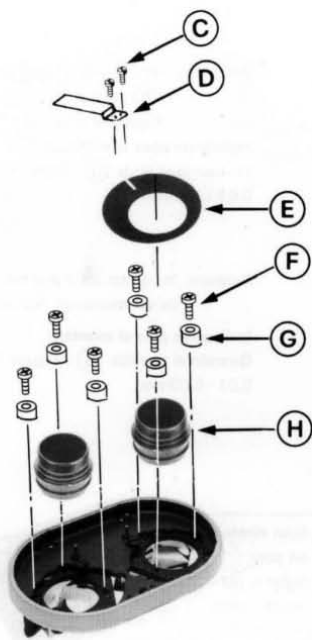Assembly instruction:

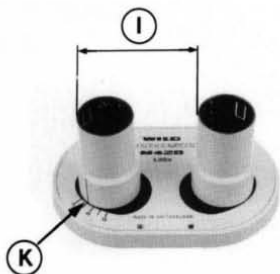
- Separating metal sheet (D) must be on top of cover disk (E).
- Measure above the tubes of the eyepieces the distance (I) and set the eye base index (K) to the measured distance by rotating the tube of the eyepiece.

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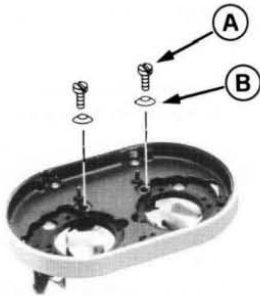
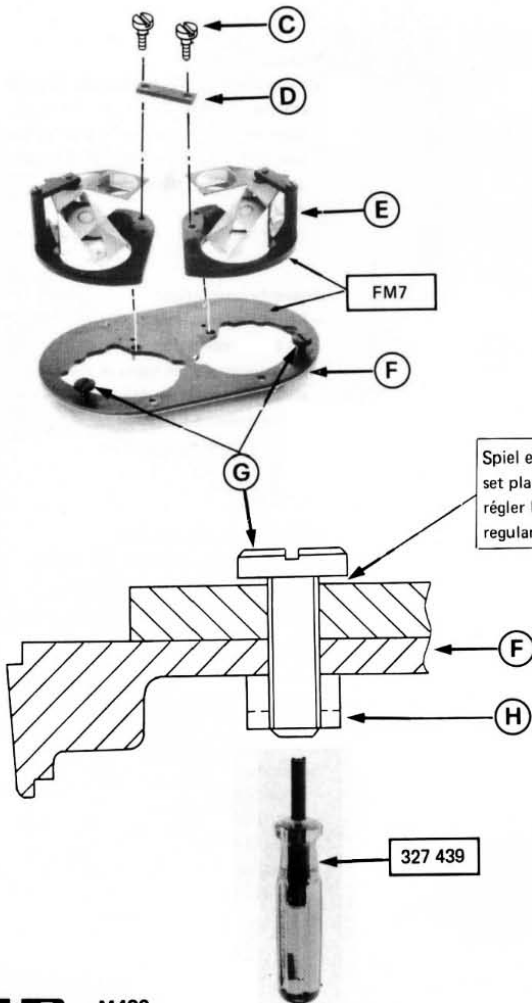
Caution: do not take the two porro prisms (E) out of their mount. Factory adjustment

Assembly instruction:

Set backlash to 0,01 to 0,03mm by rotating screw (G) .

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

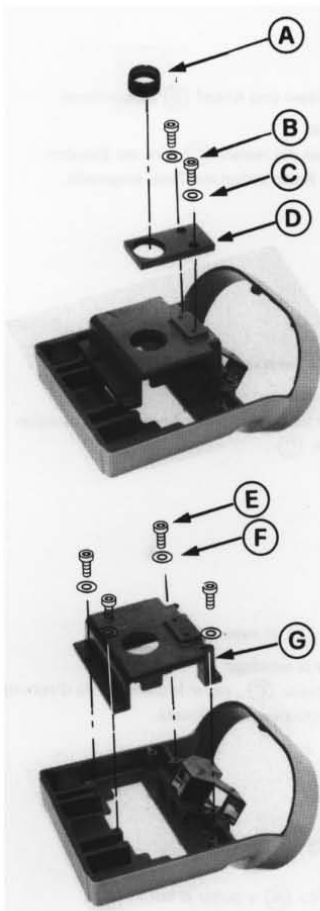
For adjusting the tubus lens (A) , the prisms (G)
and (I) see adjustment instructions

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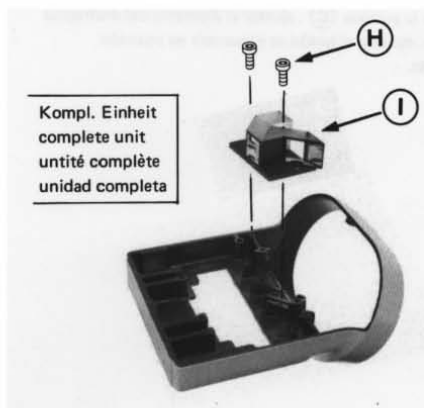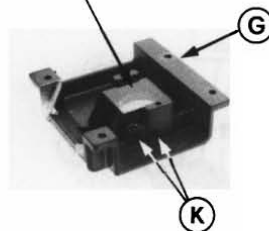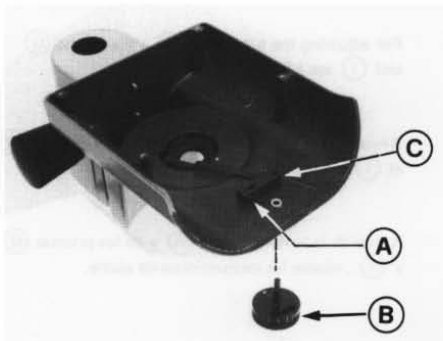
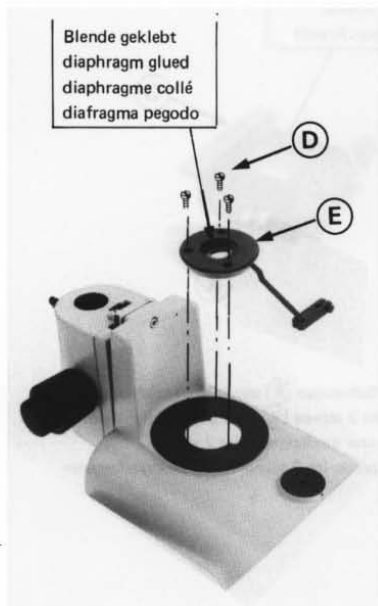
Loosen screw (A) and remove knob (B) .

Assembly instruction:

Adjust diameter of diaphragm to 7mm in latch position by rotating lever (C) .

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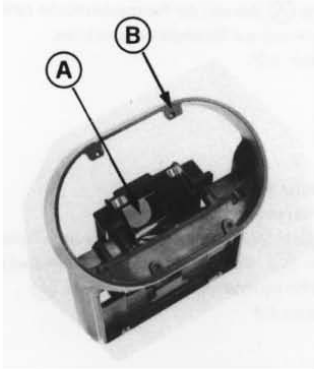
Adjusting the beam splitter prism

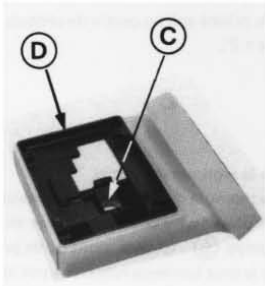
Condition:

- The distance between surface (A) of prism and the fitting surface (B) must be 73mm
- The surface (A) of prism must be parallel to the fitting surface (B)
- The surface (C) of prism must be parallel to the fitting surface (D)

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
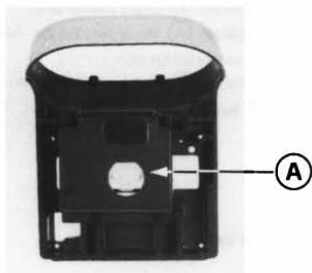
1. Adjusting the prism mount

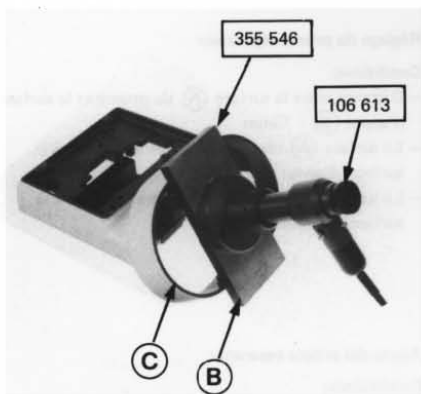- Fit prism mount (A) symmetrical
 - Place plate (B) with autocollimator onto fitting surface (C). Center the reflected illuminated cross on reticle by rotating the prism mount (A).
- Tolerance $\pm 2'$

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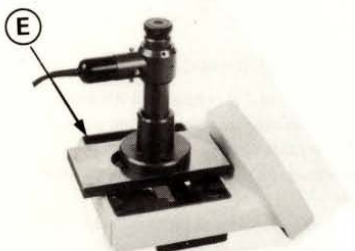
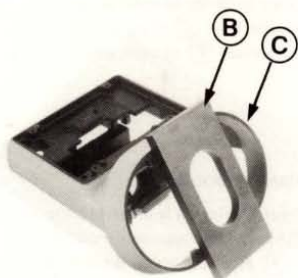
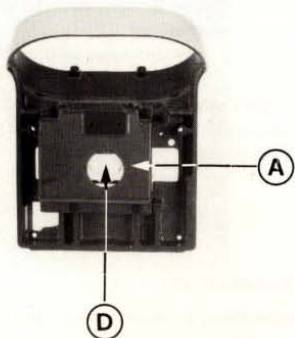
2. Adjusting the prism

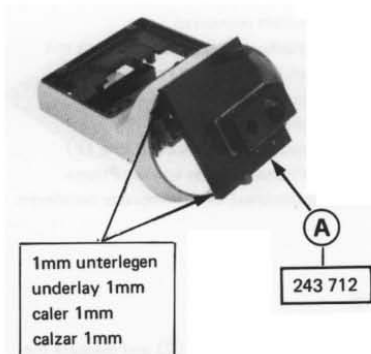
- Fit prism mount (A) symmetrical
- Place plate (B) onto the fitting surface (C) and measure the distance with the caliper. Set correct distance by shifting prism (D)
Distance = 73mm + width of plate
- Place plate (B) with autocollimator onto fitting surface (E). Center the reflected illuminated cross on reticle by tilting prism (D).
Tolerance: $\pm 5'$.

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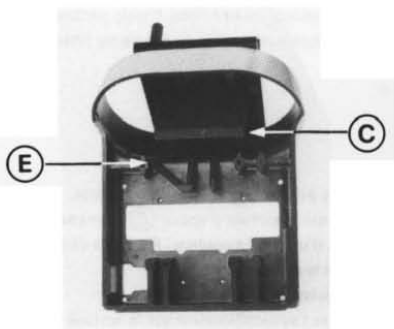
ADJUSTING THE BEAM SPLITTER PRISM

Caution: To ensure fitting the prism in the correct distance, 1mm must be underlayed between Jig (A) and optic carrier.


- Fit jig (A) with 1mm underlay to optic carrier
- Press prism (B) against edge (C) of jig and prism mount (D) against surface (E) of optic carrier

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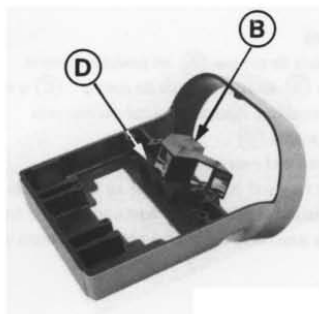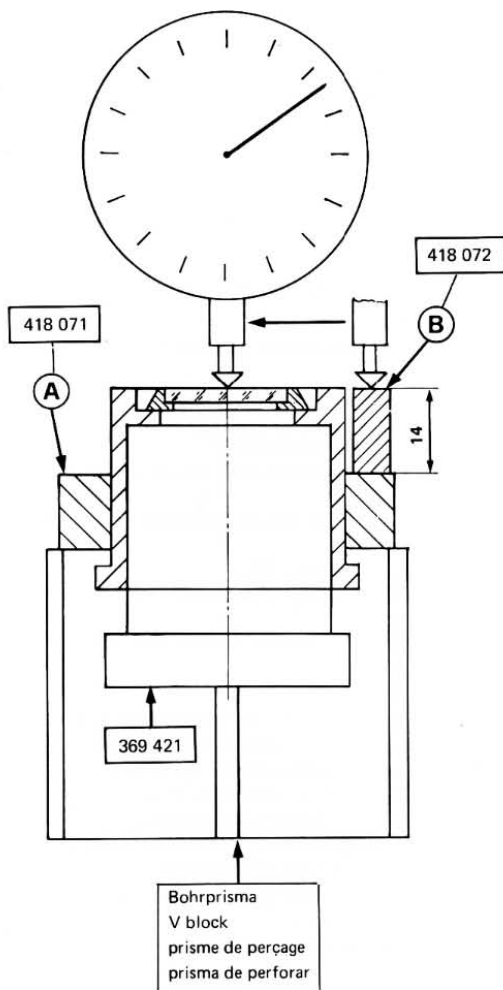
ADJUST AUXILIARY EYEPIECES

- Stick spacer ring (A) onto auxiliary eyepiece
- Set with distance gage (B) and dial gage distance of $14 \pm 0,01$ mm between reticle and reference surface
- Lock spacer ring (A).

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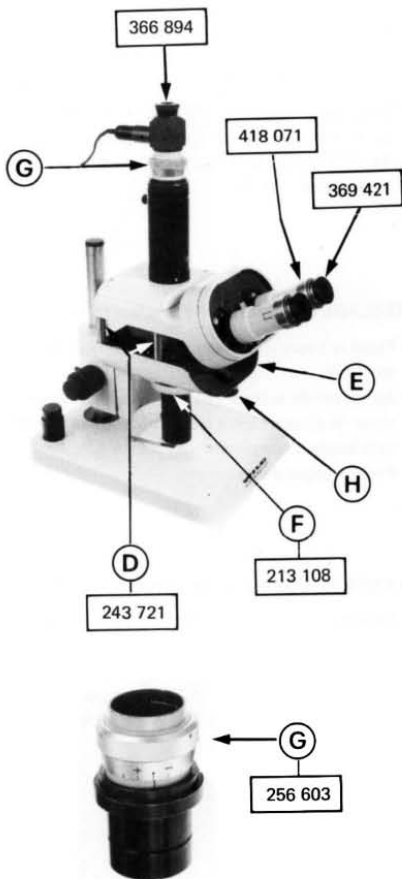
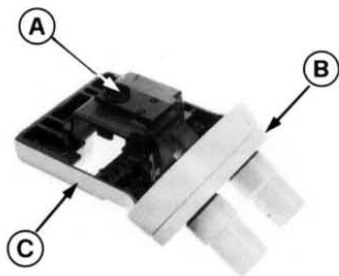
ADJUSTING TUBES LENS AND BINOCULAR TUBE

Preparation:

- Fit tube lens (A) and binocular tube symmetrical to optical
carrier (C).
- Fit optical carrier (C), separated by supports (D) to lower
part (E).
- Press mirror (F) against ring dovetail
- Insert auxiliary eyepieces with spacer ring (see pg 11) into
eyepiece tubes
- Insert adjusting tube (G) with autocollimation eyepiece into
carrier tube
- Completely open diaphragm (H).

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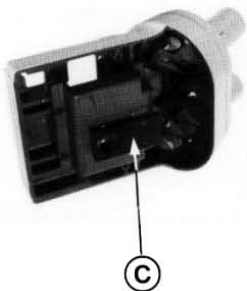
ADJUSTING TUBE LENS

1. Centering

The illuminated cross reflected from mirror (A) must be centered relative to the reticle of autocollimation eyepiece (B).

Tolerance: Small circle of autocollimation eyepiece.

Correction: Shift carrier of tube lens (C).

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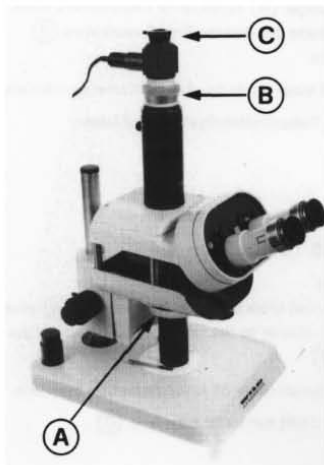
2. Height of image

The illuminated cross reflected from mirror (A) must appear sharp in reticle plane of the autocollimation eyepiece (C) when diopter setting of adjusting tube (B) is on zero.

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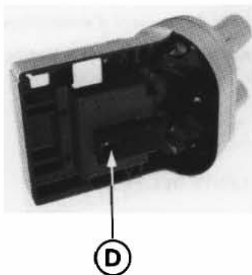
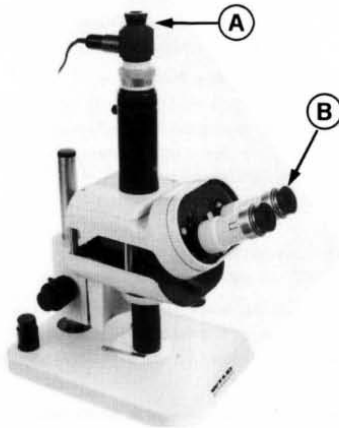
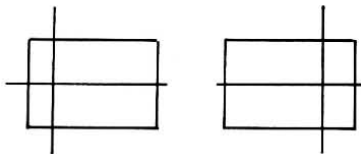
ADJUSTING BINOCULAR TUBE

1. Binocular adjustment

- The illuminated cross of the autocollimation eyepiece (A) must be within the tolerance field of the auxiliary eyepiece (B) for both eyepiece tubes.
- The difference in height from the left to the right tube may not exceed the half of the tolerance field
- These tolerances are valid for the whole range of the eyebase.

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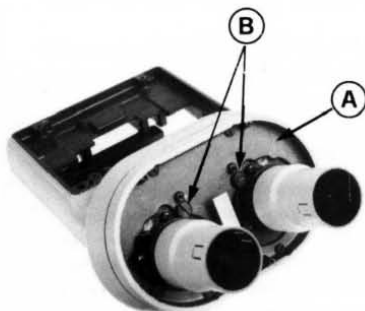
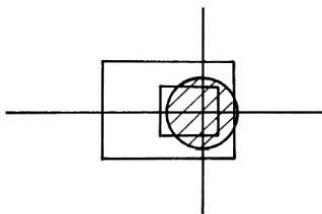
Correction of common centering errors

Shift front plate (A) parallel

Correction of binocular errors

Vertical error: rotate front plate (A)

Lateral error: loosen screw (B) and shift tube together with rotating disk.

Tolerance of binocular tube relative to photo tube:

In the right tube the reticle of the auxiliary eyepiece must be within the small circle of the autocollimation eyepiece. The middle of the photo tube must always be in the range between the two eyepiece tubes.

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2. Length of tubes

The difference between the photo tube and

- the right eyepiece tube may not exceed 0,1mm
 - the left eyepiece tube may not exceed 0,2mm
 - focus the reticle of the auxiliary eyepiece into the plane of the autocollimation eyepiece by rotating the diopter setting at the adjusting tube (A).
- Read error. 1 Interval = 0,6mm

Correction: Machine off the eyepiece tube at adjusting surface (B) on turning lathe.

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Korrektur: Abstimmung (D) abdrehen.

2. Hight of Image

The illuminated cross reflected from mirror (A) must be focused in the plane of the reticle of the autocollimation eyepiece (C) when the diopter setting at the adjusting tube (B) is on zero.

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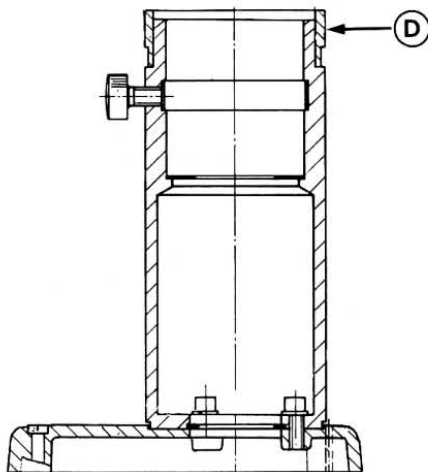
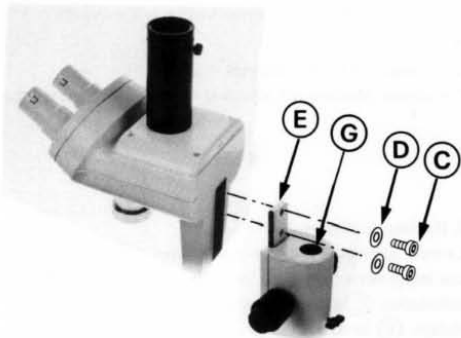
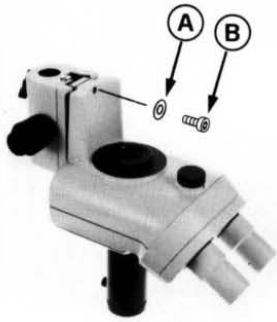
Assembly instruction:

Position hole (G) within $\pm 5'$ relative to fitting surface (H) by rotating housing (F).

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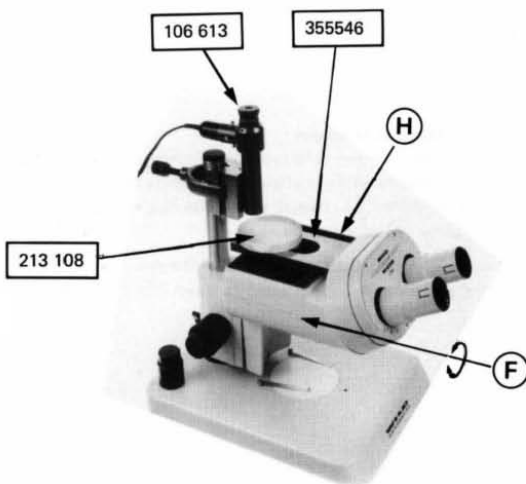
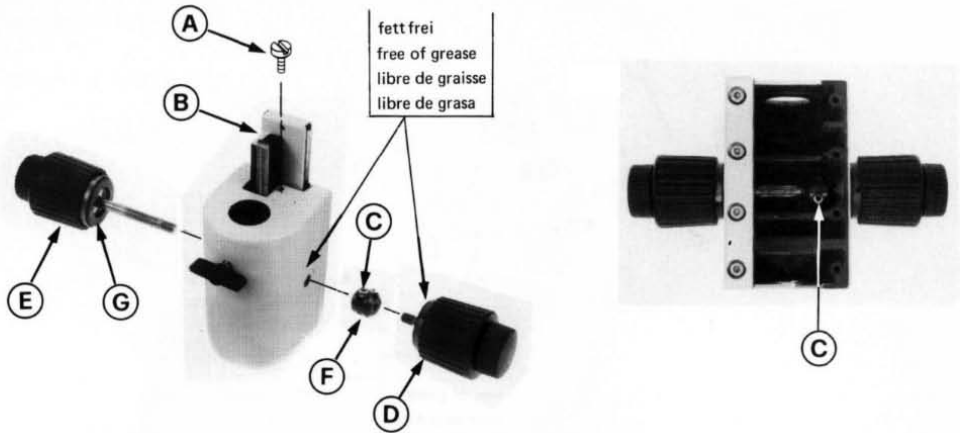
Version coarse-fine movement two-sided.

Remove screw (A) . Shift carriage (B) upwards, slacken 2 screws (C) and unscrew coarse-and fine movements (D) and (E) .

Assembly instructions:

- Clamping ring (F) must be mounted symmetrical on both spindles.
- By tightening or loosening the two coarse movements knobs adjust the braking torque of the pinion spindle with no load on the carriage to 30-40 Ncm.
- Brake disks (G) and contact surface on housing must be free of grease.

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Pinion spindle (M) is screwed into cage (K) and cemented with Araldite.

Assembly instruction:

By turning threaded ring (G) adjust braking torque to 38-42 Ncm.

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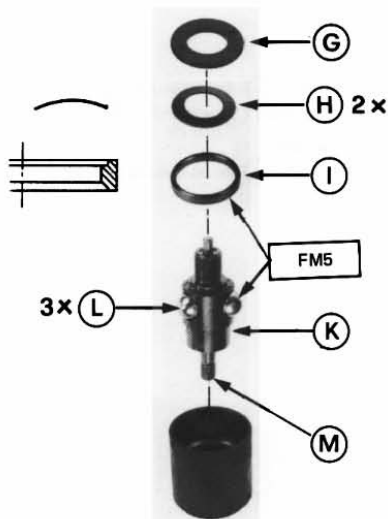
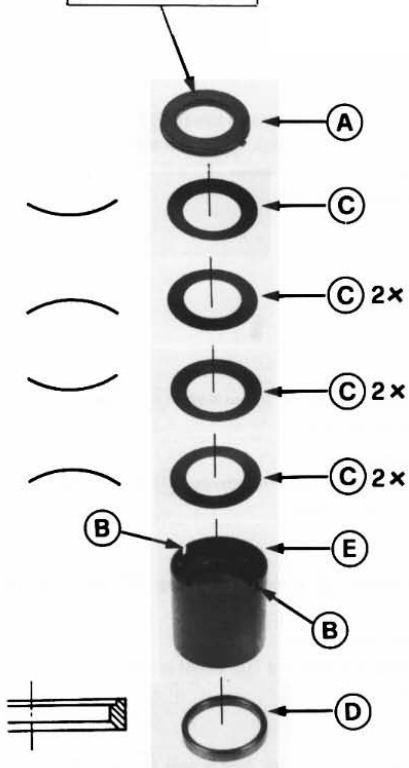
Assembly instruction: Secure brake disk (A) in the
 slots (B) of sleeve (E) .

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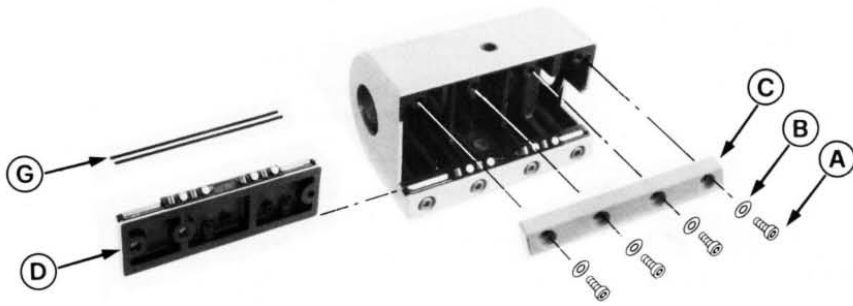
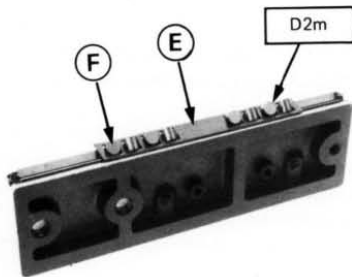
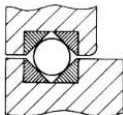
Lay housing on its side and remove screws (A), washers (B), guide bar (C), carriage (D), cages (E), rollers (F) and triangular bars (G).

Assembly instructions:

Mount rollers (F) and triangular bars (G) in correct position, see picture and drawing. Assembly of carriage see page 24

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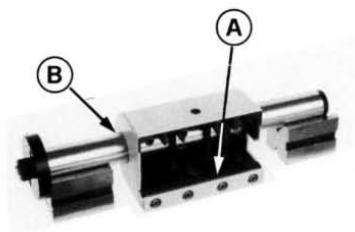
Assembly instruction:

The ball bearing track (A) must be aligned parallel to the column bore-hole (B) within 0.08mm

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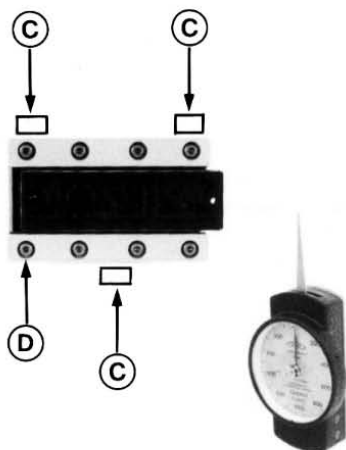
**Assembly instruction for ball track:**

The carriage must move with $5 \pm 1,5$ N, when the pinion-axis is not engaged. Mount drive housing and clamp in vise at 3 points (C) until the carriage moves with approx. 10N. Tighten screws (D).

Assembly instruction for roller track:

Mount drive housing and clamp in vise at 3 points (C) until the carriage moves with approx. 0,8-1N. Tighten screws (D).

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Assembly instruction:

- Tighten the 2 screws (B) and release again by 1/8 of a turn.
- Tighten 2 screws (A) until rack (D) rests on the screws (E).
- Set clearance between rack and pinion to 0,02 - 0,05mm by means of 2 screws (A) and (E).

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Attention:

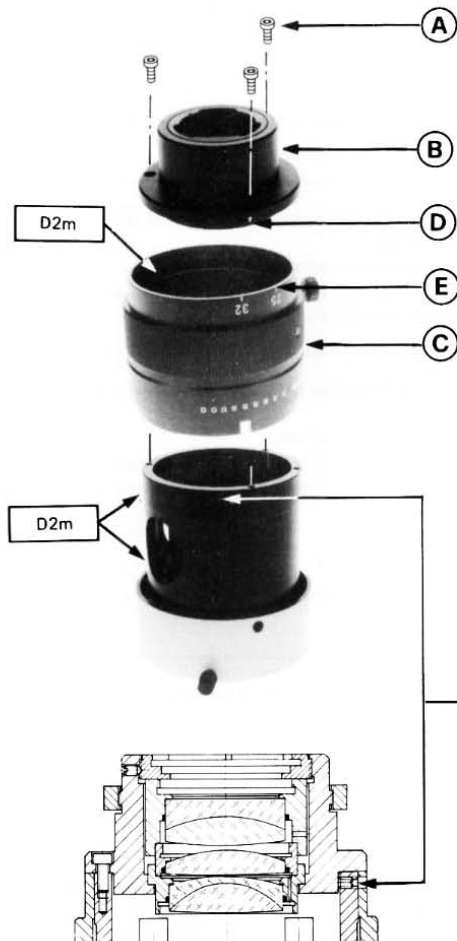
The only work which can be carried out on the objective by Service Departments is cleaning and regreasing. This is to ensure that the high quality of the objective is retained. No spare parts that would affect the adjustment will be made available.

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Attention:

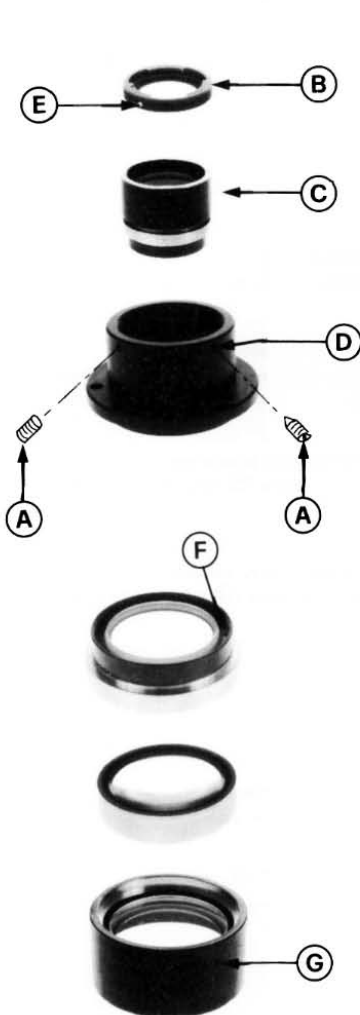
Separate the main objective only in emergency cases!

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Remove 2 screws (A) , ring (B) and objective (C) .

Mounting tip:

Tighten ring (B) until hole (D) is aligned to threaded
hole (E) .

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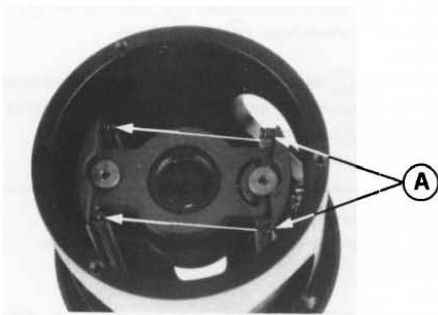
Separate objective unit with compressed air through holes
(F) and (G) .

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Place objective on the table and unhook 4 springs (A). Remove springs (A), slide (B) and (C).

Mounting tip:

Slide (C) small objective, slide (B) large objective.

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
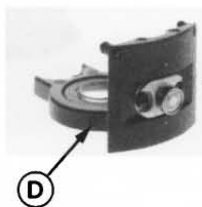
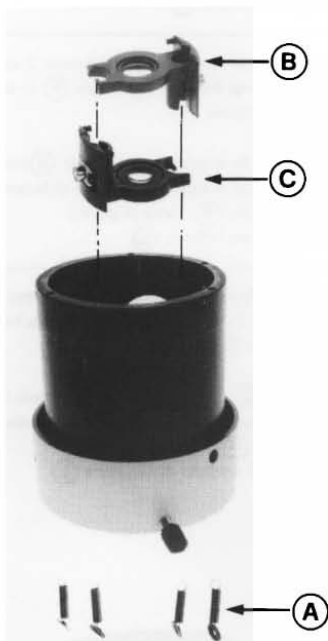
Attention: Do not dismantle optic slide!
Factory adjustment.
Do not touch adjusting screw (D)
Factory adjustment!

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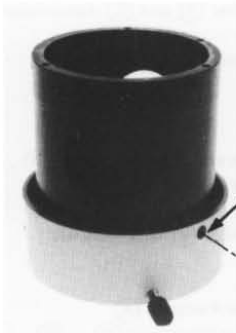Remove through hole (A) screw (B). Remove 2 screws (C), 2 springs (D), 2 brake pins (E), 3 screws (F), carrier tube (G) and camsleeve (H).

Mounting tip:

Polished surface (I) of brake pins (E) must face camsleeve (H). Orientate carrier tube (G) and tube holder (K) after slot (L) and hole (A). See photo.

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